


MESTRA®

LAVADORA POLIMERIZADORA E2
E2 WASHING AND POLYMERING MACHINE
LAVEUSE POLYMERISEUSE E2
R-080060

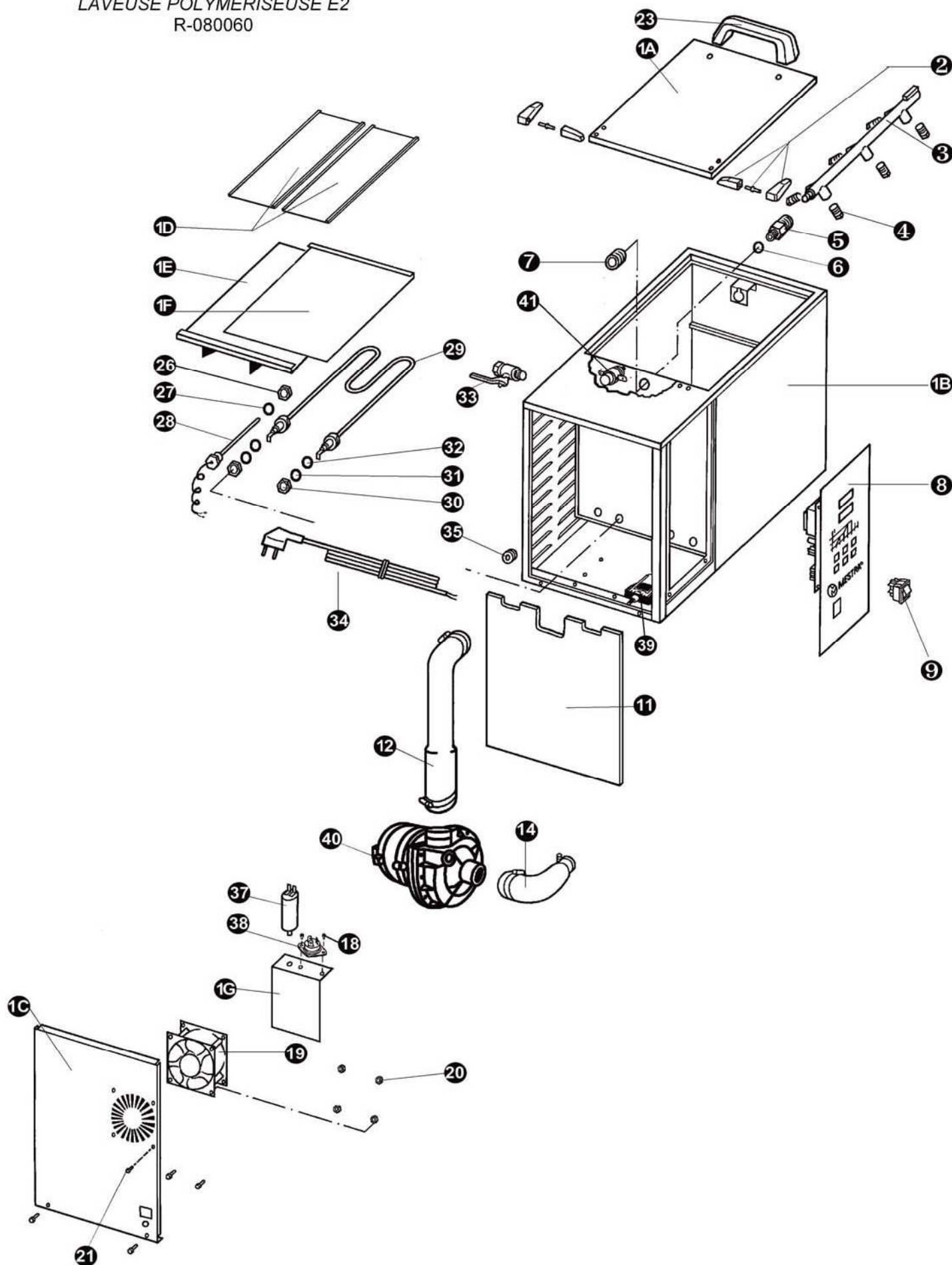

MESTRA®

LAVADORA POLIMERIZADORA E2
E2 WASHING AND POLYMERING MACHINE
LAVEUSE POLYMERISEUSE E2
R-080060 / R-080060B

Nº	Ref.	DESCRIPCIÓN	DESCRIPTION
1-A	080060-01A	Carcasa tapa superior	Housing upper lid
1-B	080060-01B	Carcasa base	Housing base
1-C	080060-01C	Carcasa lateral	Housing side
1-D	080060-01D	Bandeja estrecha	Narrow tray
1-E	080060-01E	Bandeja inferior	Lower tray
1-F	080060-01F	Bandeja ancha	Wide tray
1-G	080060-01G	Soporte de carcasa	Tray support
2	080060-02	Bisagras juego	Hinges kit
	080060-02A	Bisagras derecha	Right hinge
	080060-02B	Bisagras izquierda	Left hinge
3	080060-03	Árbol completo	Complete tree
4	080060-04	Boquilla de árbol	Tree nozzle
5	080060-05	Acoplamiento rápido	Fast coupling
6	080060-06	Arandela de cobre	Copper washer
7	080060-07	Tuerca acoplamiento manguera	Hose connecting nut
8	080060-08	Control	Control panel
9	080400-10	Interruptor 230V	Switch 230V
11	080060-11	Placa aislante	Ceramics
12	080060-12	Manguera con abrazaderas (larga)	Hose with clamp (long)
14	080060-14	Manguera con abrazaderas (corta)	Hose with clamp (short)
18	080077-10	Tornillos triac	Triac nut
19	080060-19	Ventilador	Fan
20	080490-37	Tuerca del ventilador	Fan nut
21	080075-08	Tornillo ventilador	Fan screw
22	080077-10	Tornillo de tapa de registro	Ispection cover screw
23	080400-12	Asa	Grip
26	030420-22	Tuerca de termopar	Termocouple nut
27	030420-14	Arandela de termopar	Termocouple Washer
28	080400-23	Termopar	Termocouple
29	080060-29	Resistencia	Resistor
30	080400-27	Tuerca de resistencia	Resistor nut
31	080400-26	Arandela de cobre de resistencia	Resistor copper washer
32	080400-25	Junta tórica de resistencia	Resistor O-ring
33	080400-19	Grifo de desagüe	Drain cock
34	000212	Cable 230V	Cable 230V
	000261	Cable 230V (clavija americana)	Cable 230V (American plug)
35	080070-12	Pasa cables	Cable bushing
37	080060-37	Condensador	Capacitor
38	080060-24	Triac	Triac
39	080060-41	Disyuntor	Circuit breaker
40	080060-36	Motobomba (50 Hz)	Motor pump (50 Hz)
	080060-61	Motobomba (60 Hz)	Motor pump (60 Hz)
41	080400-29	Contactador	Thermal relay