


**MESTRA®**  
POLIMERIZADORA M-18  
POLYMERIZING M-18  
POLYMERISEUSE M-18  
R-080402

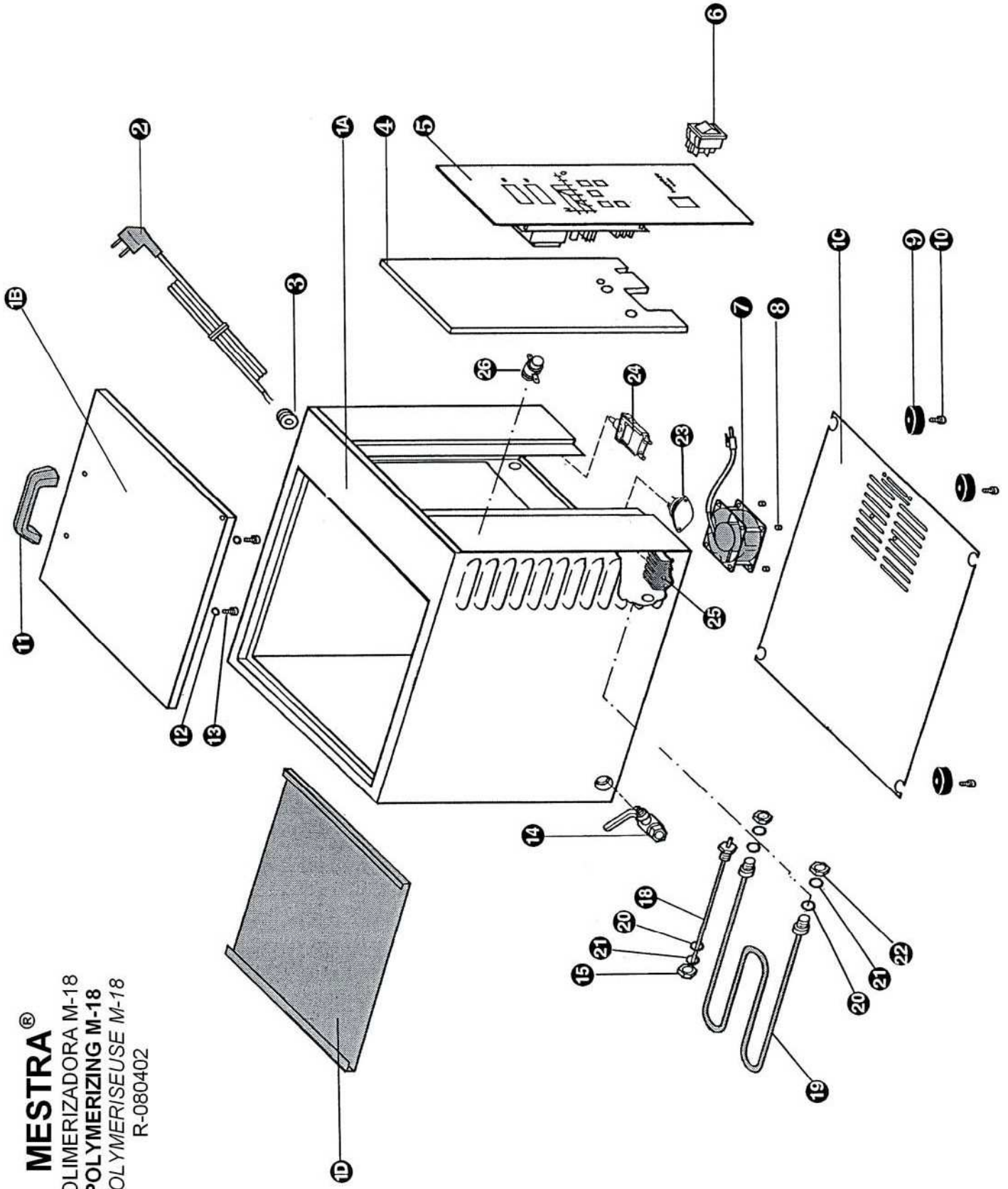

# MESTRA®

POLIMERIZADORA M-18

**POLYMERIZING M-18**

*POLYMERISEUSE M-18*

R-080402

Nº	Ref.	CONCEPTO	DESCRIPTION
1A	080402-01	Caja de acero	Stell box
1B	080402-08	Tapa superior	Upper lid
1C	080402-09	Tapa inferior	Lower lid
1D	080402-10	Bandeja interior	Inner salver
2	000212	Cable	Cable
	000214	Cable (América)	Cable (America)
3	080010-05	Pasa cables	Whire bushing
4	000302	Cerámica	Ceramic
5	080400-06	Control	Control panel
6	080400-10	Interruptor	Switch
7	080400-08	Ventilador	Fan
8	080400-09	Casquillo	Bushing
9	080080-26	Patas	Legs
10	080251-09	Tornillos de patas	Legs screws
11	080400-12	Asa	Handle
12	080460-13	Arandela de asa	Handle washer
13	080460-65	Tornillo de asa	Handle screw
14	080400-19	Grifo de desagüe	Drainpipe tap
15	030420-22	Tuerca termopar	Thermocouple nut
18	080400-23	Termopar	Thermocouple
19	080402-04	Resistencias	Resistor
20	080400-25	Junta tórica resistencia	Resistor O-ring
21	030420-14	Arandela resistencia	Resistor washer
22	080400-27	Tuerca de resistencia	Resistor nut
23	080060-24	Triac	Triac
24	080060-41	Disyuntor	Circuit breaker
25	080402-06	Disipador	Heat sink
26	080400-29	Contactador	Contactador