


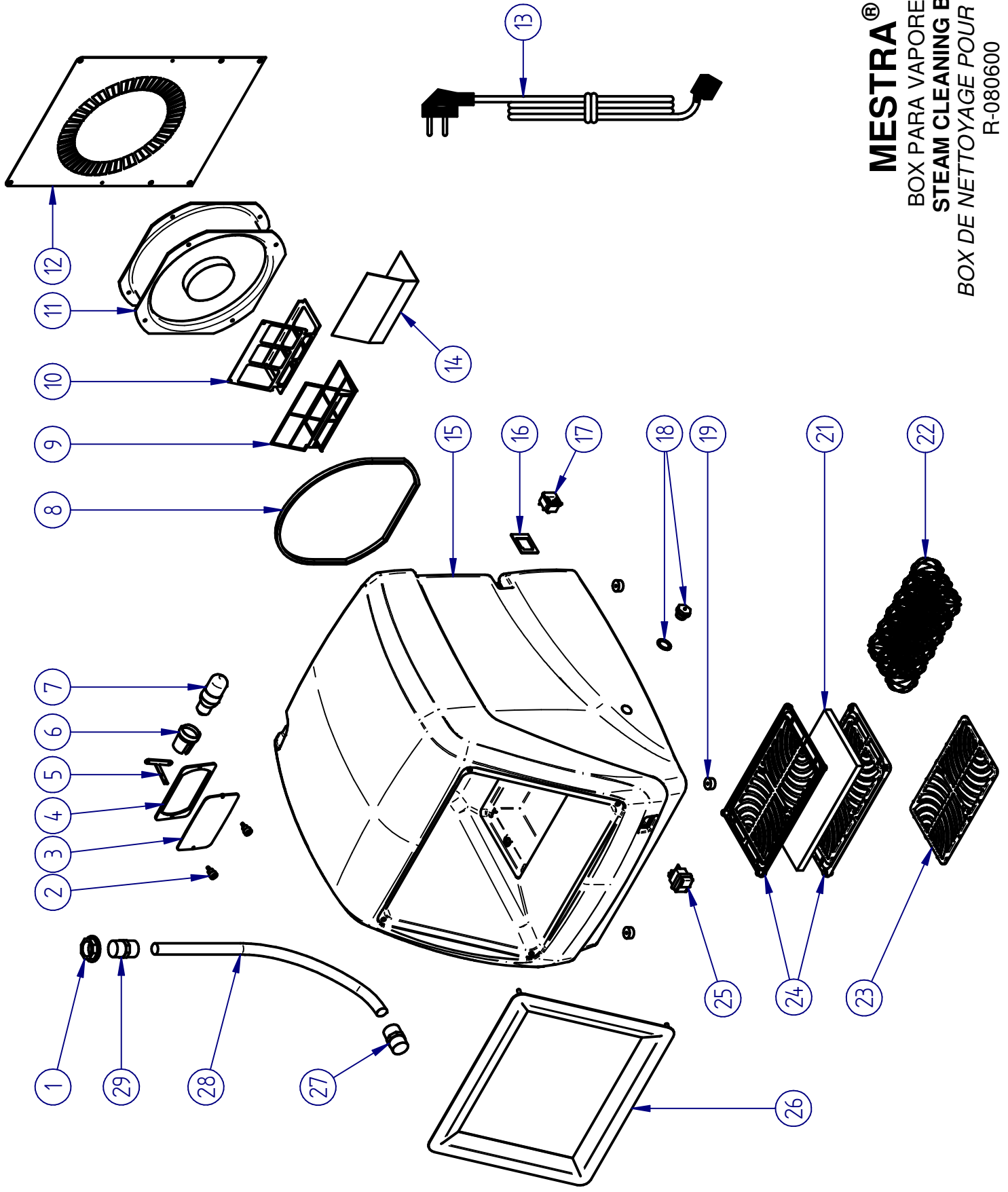

MESTRA®
 BOX PARA VAPORETO
 STEAM CLEANING BOX
 BOX DE NETTOYAGE POUR LA VAPEUR
 R-080600

MESTRA®
 BOX PARA VAPORETO
STEAM CLEANING BOX
BOX DE NETTOYAGE POUR LA VAPEUR
 R-080600

Nº	Ref.	CONCEPTO	DESCRIPTION
1	080086-06	Tuerca de manguito	Sleeve nut
2	080600-28	Pomo pantalla	Shield knob
3	080600-02	Pantalla lámpara	Lamp shield
4	080600-07	Junta pantalla	Shiels gasket
5	080600-16	Soporte portalámparas	Lamp holder support
6	080600-13	Portalámparas	Lamp holder
7	080600-05	Lámpara 230V	Lamp 230V
	080600-05B	Lámpara 115V	Lamp 115V
8	080600-26	Junta de ventilador	Fan gasket
9	080600-19	Marco filtro ventilador	Fan filter frame
10	080600-20	Empujador filtro ventilador	Fan filter pusher
11	080600-03	Ventilador 230V	Fan 230V
	080600-03B	Ventilador 115V	Fan 115V
12	080600-18	Chapa trasera	Back sheet
13	000250	Cable 230V	Cable 230V
	000252	Cable 115V	Cable 115V
14	080600-21	Filtro de espuma	Foam filter
15	080600-08	Carcasa	Housing
16	080600-17	Soporte base porta fusibles	Fuse holder base support
17	080490-06	Base porta fusibles	Fuse holder base
18	080600-01	Visor	Visor
19	080490-31	Patas	Legs
21	080600-22	Filtro estropajo	Cloth filter
22	080600-12	Conjunto filtros metálicos	Metal filter kit
23	080600-40	Rejilla inferior cámara	Chamber lower grid
24	080600-39	Rejillas superiores de cámara	Chamber top grids
25	080400-07	Interruptor	Switch
26	080600-09	Marco frontal	Front frame
27	080086-07	Manguito inferior de manguera	Lower hose sleeve
28	080600-04	Manguera de desagüe	Hose
29	080210-02	Manguito superior manguera	Top hose sleeve