


Spare parts 18030500 Version A



Reparaturzeichnung • Repair services drawings • Eclatés pour les services de réparation • Esplosi per servizi d'assistenza
 Diseños para servicios de reparaciones • Загрузка чертежей Для ремонтных мастерских

1803 xxxx MT plus

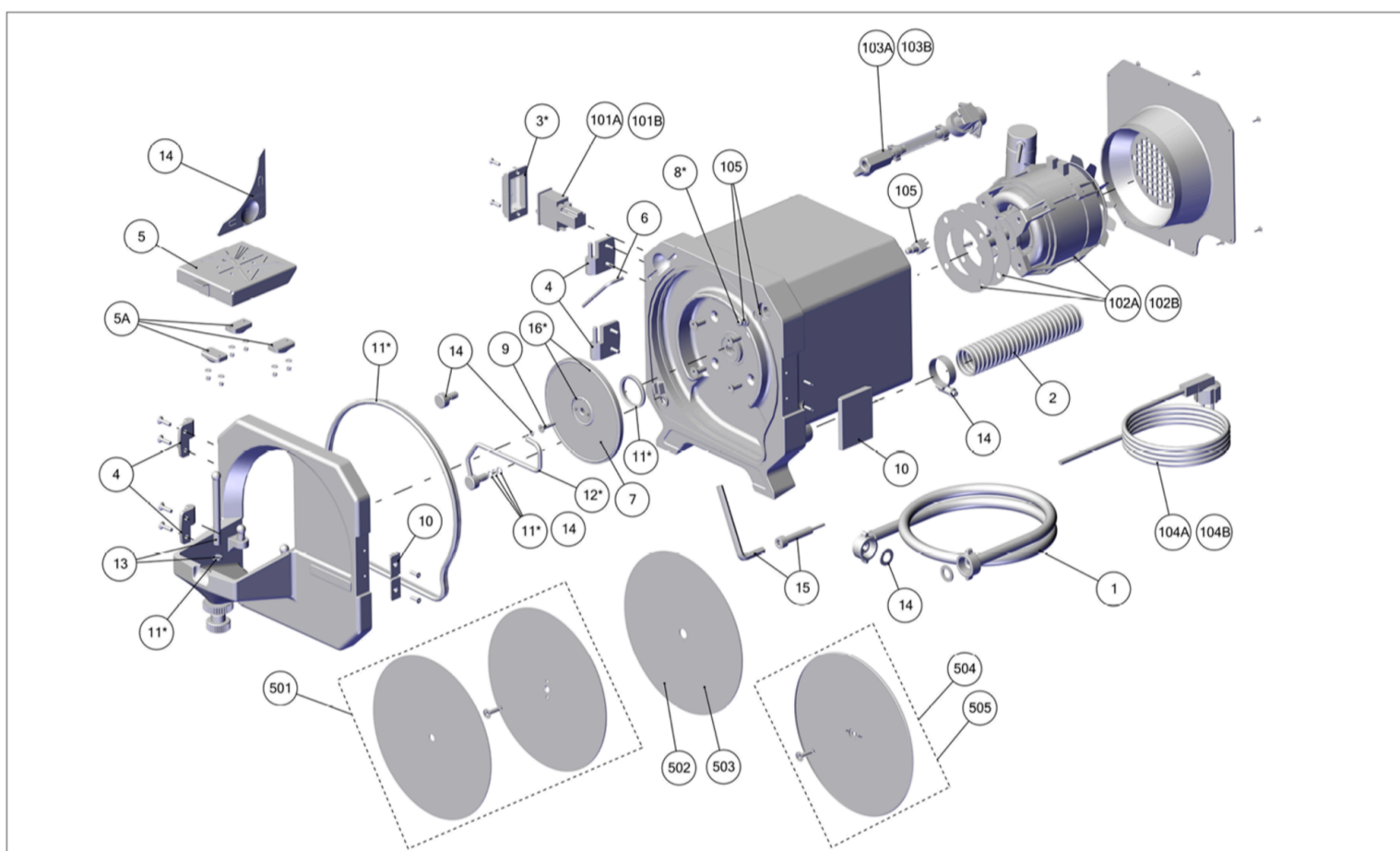

21-9874 18012018 / A

Position	Product name	Item number
1	Admission hose compl. Trimmer	20976
2	Outlet hose 41x35mm grey 1,5m long	900150449 *
3	Protection cap on/off switch MTplus compl.	900035788 *
4	MTplus Hinges complete	900035708
5	MTplus Trimmer table complete	900035717
5A	Ballsocket Trimmer table	900021357
6	Hexagon-key SW 4	900033137
7	Flange MTplus complete up to stand B (O-ring)	900033454



Position	Product name	Item number
8	Protection cap door switch compl.	21234 *
9	Screw for trimming disc MT2/TT2/MTplus/MT3/MT3 pro	900035237
10	Door locker MTplus complete	900033150
11	Set of joints MTplus	900035711 *
12	Spray tube MTplus compl.	900035716 *
13	Adjusting screw MTplus complete	900036628
14	Small part set MT plus	900021336
15	Set of tools for MTplus	900021337
16	Set of O-rings for flange MTplus/MT2/TT2/MT3/MT3 pro	900033450 *
101A	On/Off-switch MTplus 230 V/240 V_12 A	900038698
101B	On/Off-switch MTplus 100V / 120V	900035736
102A	Driving motor MTplus 230V compl.	900035719
103A	Water admission MTplus 230V / 240V compl.	918039998
103B	Water admission MTplus 100V / 120V compl.	918039997
104A	Mains cable /Harness MTplus 230V	900035720
104B	Mains cable /Harness MTplus 120V	900035744
105	Door switch Trimmer	900031881
501	Klettfix MT plus Dia. 234mm	18031000
502	Klettfix MTplus, MT3 and MT3 pro, refill pack grain 80	18031100
503	Klettfix MTplus, MT3 and MT3 pro, refill pack grain 120	18031200
504	Marathon MT plus Dia. 234mm	18032000
505	Infinity MT plus Dia. 234mm	18033000

* Wear and tear parts, excluded from the guarantee.



