


Spare parts 29342000 Version A



Reparaturzeichnung • Repair services drawings • Eclatés pour les services de réparation • Esplosi per servizi d'assistenza
Diseños para servicios de reparaciones • Загрузка чертежей для ремонтных мастерских

2934 2X00 / 2934 3X00 SILENT compactCAM

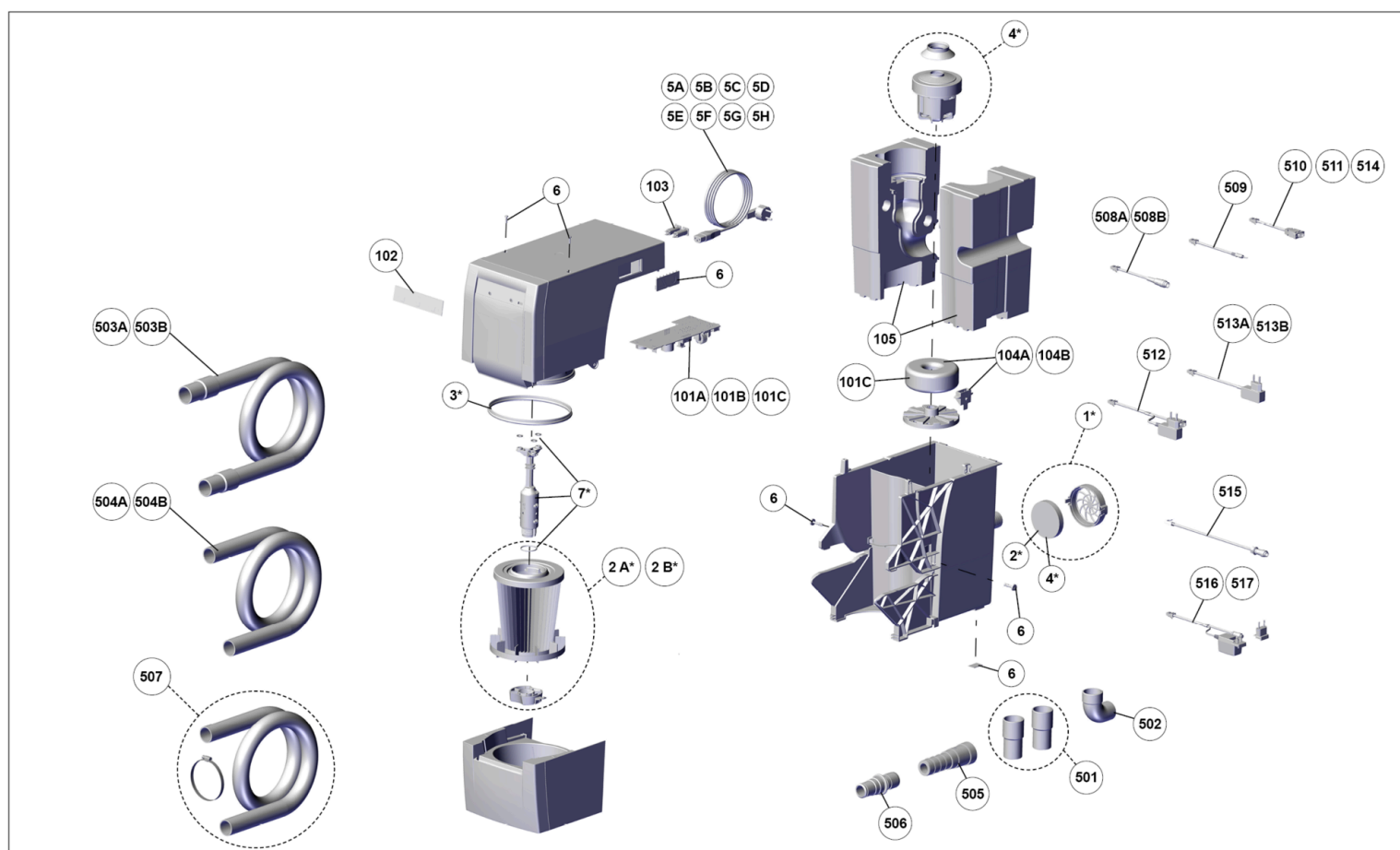

21-6831 03092024 / A

Position	Product name	Item number
1	Exhaust air filter kit	900037585 *
2A	Fine filter SILENT compact / SILENT compactCAM	900021635 *
3	Bellow seal	900037473 *
4	Turbine SILENT compact	900037226 *
5A	Mains cable 230 V with low-power plug	900032750
5B	Mains cable 120V with low-power plug	900035570
5C	Unit connector cable 230V UK	900035938
	Power cord Argentina with cable-mount connector, straight	900032769



Position	Product name	Item number
5E	Unit connector cable 230V CH	900036070
5F	Power cord Brazil with cable-mount connector, straight	900032767
5G	Power cord SA 10A with cable-mount connector, straight	900032765
5H	Power cord Italy with cable-mount connector, straight	900032768
6	Kit of small parts	900037238
7	Vibration motor for cleaning	929349998 *
101A	Printed board circuit SILENT compactCAM, 230V / 220V	900033514
101B	Printed board circuit SILENT compactCAM, 120V / 100V	900033515
101C	Printed board circuit SILENT compactCAM, 220V (only stand A)	900037514
102	Operating foil SILENT compactCAM	900037462
103	Appliance protective switch 8A	900037389
104A	Toroidal core transformer 230VAC / 120VAC	900037224
104B	Toroidal core transformer 220 VAC / 100 VAC	900037682
105	Foam inlay complete	900021590
501	End bushing set, 2 pieces	29210003
502	90° angled connector Silent	29340007
503A	Suction Hose	900034240
503B	Suction hose, antistatic, 3 m, incl. 2 end mufflers	900034826
504A	Suction hose inside diameter 38 mm, 6 m	901150823
504B	Suction hose inside diameter 38 mm, 9 m	902150823
505	Adapter for hose connection	900034305
506	Universal suction hose adapter	900034430
507	External exhaust air adapter SILENT compact	29340004

Position	Product name	Item number
508A	Interface cable Type A	29340005
508B	Interface cable Type G	29340013
509	Interface cable Type B	29340006
510	Interface cable Type C	29340008
511	Interface cable Type D	29340009
512	Interface cable Type E + Nema adapter	29340010
513A	Interface cable Type F	29340011
513B	Interface cable Type F + C14 adapter	29340012
514	Interface cable Type H	29340016
515	Interface cable type I	29340017
516	Interface cable type J	29340018
517	Interface cable type M	29340021

* Wear and tear parts, excluded from the guarantee.

