


Spare parts 29590000 Version A



Ersatzteilzeichnung • Spare parts drawings • Plan des pièces de rechange • Disegno delle parti di ricambio
Plano de piezas de recambio • Чертеж запасных частей

2959-xxxx Basic quattro IS

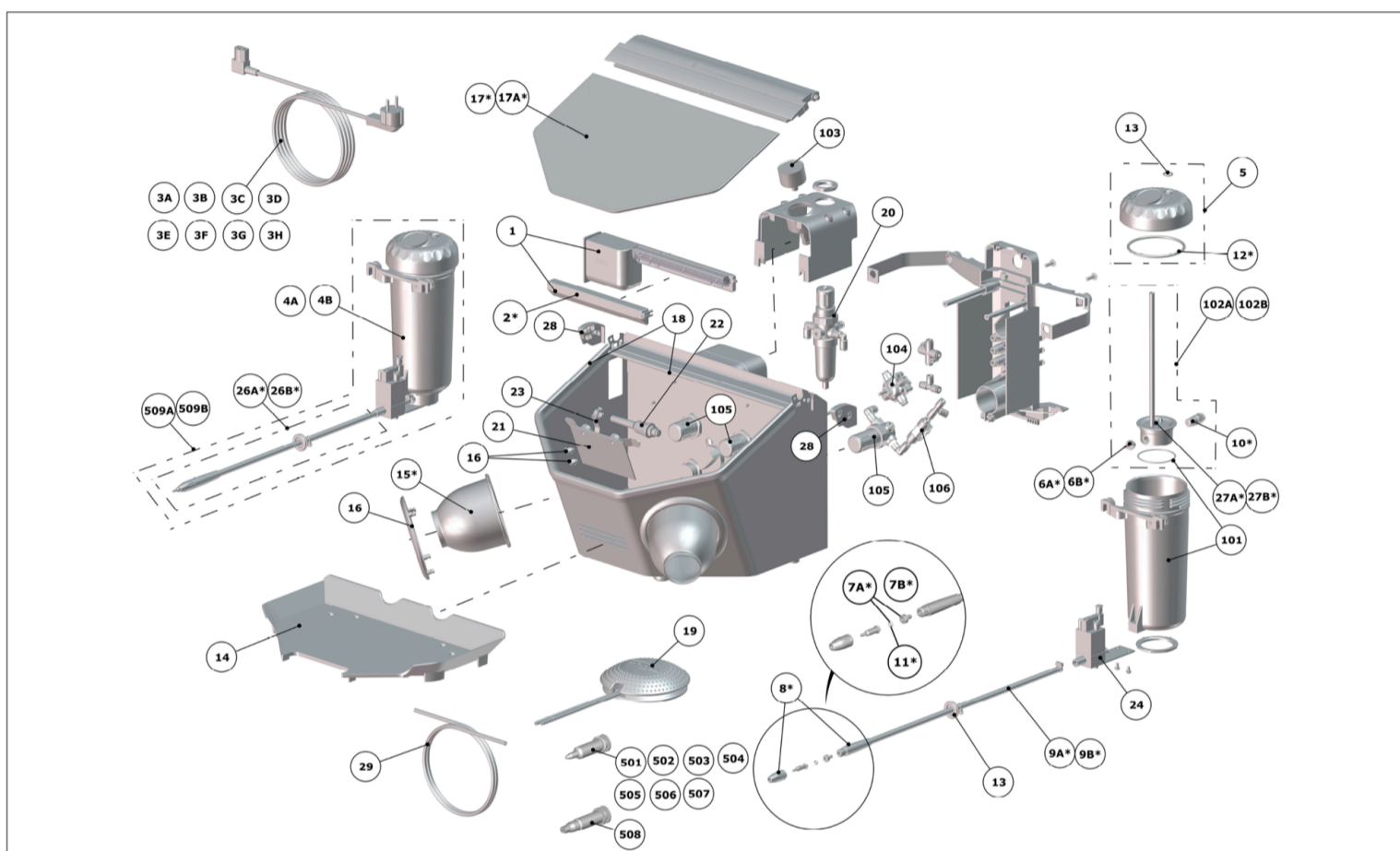

21-5644 20102025 / A

Position	Product name	Item number
1	Lamp complete	900021428
2	Cover for lamp	900036659 *
3A	Unit connector cable 230V	900036121
3B	Unit connector cable 120V	900036122
4A	Replacement tank Basic IS 25-70µm	900021455
4B	Replacement tank Basic IS 70-250 µm	900021456
5	Tank lid IS	900036398
10	Filter cartridge	900021431 *
	O-ring DIN 3771 82x4 NBR 55	900033035 *



Position	Product name	Item number
13	Hardware Basic	929470002
14	Base grid	900036330
15	Hand sleeves	900034751 *
16	Ring for hand sleeves	900036339
17	Coated glass screen master / quattro	900021435 *
17A	Uncoated glass screen master / quattro (3 pcs.)	900036311 *
18	Edge protection Basic	929470001
19	Pneumatic foot switch	900020921
20	Maintenance unit IS	900036600
21	Ventilating flap 2946	900034104
22	Air nozzle, complete	929540001
23	Hose feed fixation quattro / quattro IS	900036683
24	Valve unit IS	900036409
26A	IT blasting hose incl. handpiece 25-70 µm	929000050 *
26B	IT blasting hose incl. handpiece 70-500 µm	929000500 *
27A	Plug-in dosing nipple 25-70 µm	900036353
27B	Plug-in dosing nipple 70-250µm	900036354
28	Side panels Basic	929470003
29	PU-hose 6 x 3.9 mm, blue 5 M	900150864
101	Replacement tank case Basic	900035800
102A	Mixing chamber IS 25-70µm Basic quattro IS	900036421
102B	Mixing chamber IS 70-250µm Basic quattro IS	900036422
103	Manometer with screw fitting	21119
104	4-fold tank selector switch	900021550
105	Pressure regulator IS Basic quattro IS	900036414
106	Check valve IS	900036415

Position	Product name	Item number
501	IT Blasting nozzle 0,4 mm	900021203 *
502	IT Blasting nozzle 0,6 mm	900033213 *
503	IT Blasting nozzle 0,8 mm	900021204 *
504	IT Blasting nozzle 1,0 mm	900033211 *
505	IT Blasting nozzle 1,2 mm	900033214 *
506	IT Blasting nozzle 1,4 mm	900021205 *
507	IT Blasting nozzle 2,0 mm	900021206 *
508	IT Slotted nozzle 1,5x3,5 mm	900031739
509A	Additional tank 25-70 μ basic quattro IS 2959	29590050
509B	Additional tank 70-250 μ basic quattro IS 2959	29590250

* Wear and tear parts, excluded from the guarantee.